

How to save your distance measurement results in DV Softworx

1. drag data to window here

2. select 2-point distance

3. click save options

4. click "enter" afterwards

5. tick

6. click "save distance"

7. "measure" distance

8. back to the image, select all points for distances. It will auto save all distances in the folder defined in step 4

The screenshot shows the DV Softworx 4.1.0 interface. The 'Measure Distances' window is open, showing 'Standard Two Point' as the measurement method and 'Micrometers' as the units. The 'Distance Save Options' dialog is also open, with the 'Single Column' checkbox checked. The background image shows a grid of points with lines connecting them, representing distance measurements. The taskbar at the bottom shows the software running on a host named 'DV-Analysis'.